

HAMILTON SYRINGE TIPS

RINSING, PACKING AND SHIPPING PROCEDURES

Items Supplied

- 1 – Shipping Carton with Cover Tray
- 2 – Rinsing & Shipping Containers With Solid Yellow Lids
- 2 – Drain Lids (First Shipment Only)
- 2 – Bottles Sodium Hypochlorite (Bleach)
- 1 – FedEx Ground Return Record

PROTECTIVE GLOVES SHOULD BE WORN DURING RINSING PROCESS

1. LOAD

At the end of shift, or when used syringe tray becomes full, empty used syringes from the instrument waste tray into the Yellow Lid Rinsing & Shipping Container provided.

2. FILL

(Loaded container needs to be taken to a convenient rinsing station)
Empty ½ bottle of Sodium Hypochlorite (Bleach) into the rinsing container and fill approximately one-half full with D.I. water as indicated. (If D.I. water is not conveniently available, tap water will do.)

3. RINSE

Seal container with solid lid. Hold over the sink and invert 3 or 4 times so that all syringes are exposed to the liquid and let stand for 5-10 minutes.

4. DRAIN

Replace Solid Lid with Drain Lid and pour solution down drain. Inverting the container a second and third time will allow for additional liquid to be shed from syringes. After draining, replace Drain Lid with Solid Lid for shipping.

5. BOX

Place Rinsing & Shipping Container with decontaminated syringes into bottom of Shipping Carton. Enclose empty bleach bottles in bottom portion of box also. Place tray in box and press down to be in contact with Yellow top Shipping Containers. Place syringe racks and end clips in tray along with completed shipping form.

6. SHIP

Seal Shipping Carton with preprinted return label appearing on outside flap. Call FedEx Ground at **1-800-937-4587** for pick up.

NOTE: PLEASE RETAIN DRAIN LIDS FOR FUTURE USE.

We appreciate your cooperation in following these procedures and look forward to serving you.

